


Complete bisulfite conversion of GC-rich DNA in less than 3 hours

EZ DNA Methylation-Gold® Kits

- **Consolidated Workflow:** DNA denaturation and bisulfite conversion process in one step.
- **High Yield:** Desulfonation and purification on a spin-column allows elution of bisulfite-converted DNA in as little as 10 µl.
- Suitable for Various Downstream Applications: Bisulfite-converted DNA can be robustly PCR amplified for downstream analyses including endonuclease digestion, sequencing, microarrays, etc.


PCR-Ready Bisulfite Converted DNA

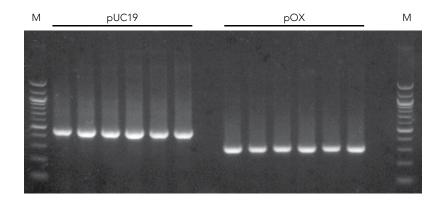

Salmon sperm DNA were spiked with fully (CpG) methylated, linearlized pUC19 and pOX plasmid (1.5% w/w) and subjected to bisulfite conversion using the EZ DNA Methylation – Gold $^\circ$ kit. Two amplicons, 453 bp and 314 bp, were successfully amplified from the bisulfite converted pUC19 and pOX plasmid DNA, repectively. The amplicons were analyzed in a 2.0% (w/v) TAE/agarose/ EtBr gel. The size marker "M" is a 100 bp ladder.

Product	Cat. No.	Size
EZ DNA Methylation-Gold®	D5030	50 rxns
EZ-96 DNA Methylation-Gold® (shallow-well)	D5032	2 x 96 rxns
EZ-96 DNA Methylation-Gold® (deep-well)	D5033	2 x 96 rxns
EZ-96 EZ DNA Methylation-Gold® Magprep	D5046	4 x 96 rxns

Streamlined Workflow

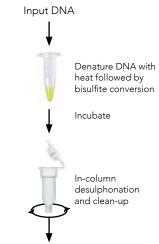

Elute Bisulfite-Converted DNA

Purified DNA is added directly to the CT-conversion reagent. Denaturation and bisulfite conversion are completed in a single-tube reaction. After desulphonation and purification in a spin-column, the purified bisulfite-converted DNA is ready for PCR.

Specifications	
Input	500 pg - 2 µg purified DNA
Conversion Efficiency	> 99%
Elution Volume	Spin Column: ≥ 10 μl 96-well: ≥ 15 μl Magprep: ≥ 25 μl
DNA Recovery	> 75%
Platform Compatibility	Swift Biosciences [™] Accel-NGS [®] Methyl-Seq DNA Library Kit
	Agilent SureSelectXT Methyl-Seq Target Enrichment System



info@zymoresearch.com



www.zymoresearch.com



Toll Free: (888) 882-9682