


Capture the Magic...

Novel Surface Chemistry for Superior cfDNA Isolation


MAGicBead™ cfDNA Isolation Kit

Novel magbead surface technology moves away from traditional silica-based DNA binding chemistry. It uses innovative DNA binding and release mechanism that is quick and efficient, free of use of alcohol.

- **Highest Yield:** Robust cfDNA isolation from up to 10mL plasma w/ unique MAGicBead™ technology
- **Compatible:** Any open automation platforms and a variety of biofluids, including plasma
- **NGS-Ready:** High-quality cfDNA ideal for Next-Gen Sequencing and other applications

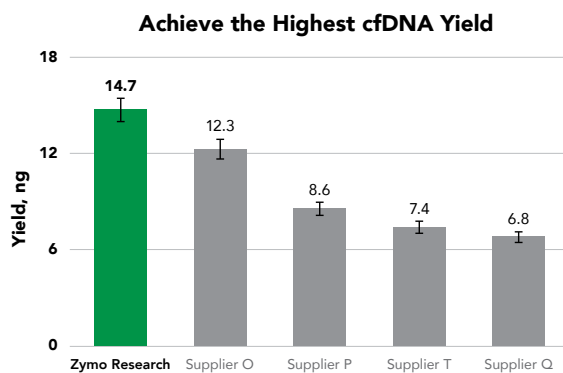

Total cfDNA yields from five different magbead-based kits. Triplicate preps of 1mL plasma each from a single donor was used. Yields were assessed using Qubit™ 1x dsDNA HS Assay Kit (Thermo-Fisher Scientific).

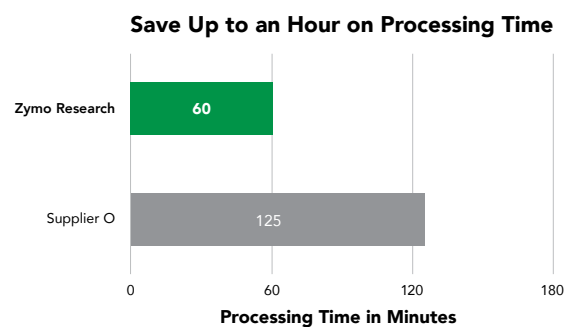


Total processing time required from the latest cfDNA isolation technologies. Plasma cfDNA was extracted from 1mL of plasma using two different magbead-based kits.



ZYMO RESEARCH


The Beauty of Science is to Make Things Simple®

Simple Workflow

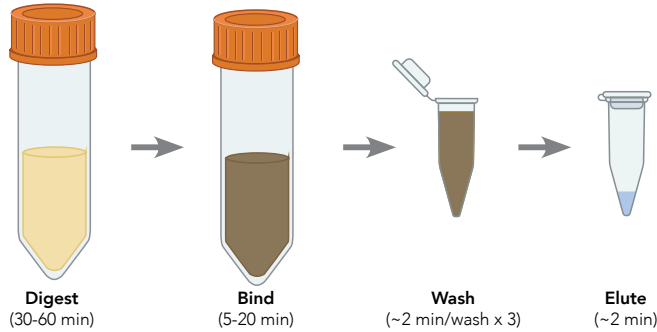

Simple, scalable, and quick protocol that is easy to adopt on any automation platform or on lab bench.


Scalable Recovery

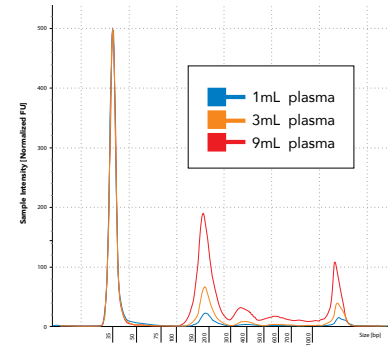

Efficiently scale your sample input from 0.2 to 10 mL

Compatibility

Sample types	Blood tube types (plasma)	Automation platforms	Incompatible
<ul style="list-style-type: none"> Plasma Serum Saliva Urine Cerebrospinal fluid Amniotic fluid Spent cell culture media 	<ul style="list-style-type: none"> K₂EDTA Streck-DNA K₃EDTA Na₂EDTA Na-Citrate Na-F / K-Oxalate 	<ul style="list-style-type: none"> KingFisher™ Flex/Apex (Thermo-Fisher) Fluent® X (Tecan) Microlab® STAR™ (Hamilton) OT-2 (Opentrons) 	<ul style="list-style-type: none"> Heparin Synovial fluid

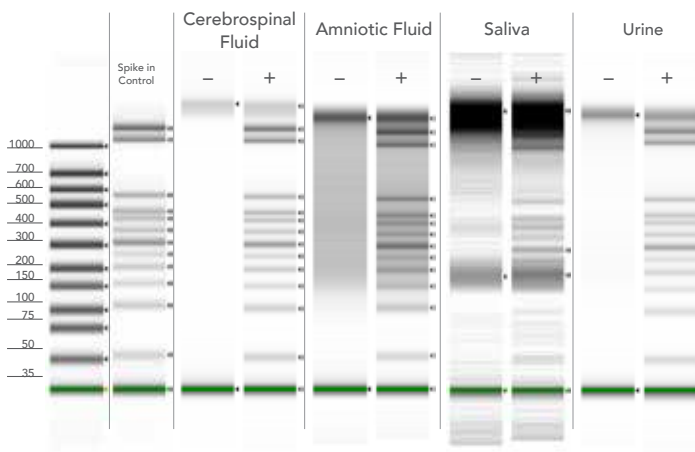

Compatible with wide variety of sample types
500 µL of cerebrospinal fluid, amniotic fluid, spun-down saliva (12,000 x g for 5 min at RT) or 3 mL of urine.

Visualized using Cell-free DNA ScreenTape Analysis (Agilent Technologies).

Experience the Magic for Free!



Product	Cat. No.	Size
MAGicBead™ cfDNA Isolation Kit	D4086	2 mL x 50 Prep.

Email automation@zymoresearch.com for assistance, including scripts and other specific inquiries tailored to your project. Require a Custom Solution? [Inquire Here](#) or email busdev@zymoresearch.com



info@zymoresearch.com



www.zymoresearch.com



Toll Free: (888) 882-9682