

Optimized Lysis Protocols

The following conditions with different mechanical lysis machines were validated with minimum bias using the ZymoBIOMICS™ Microbial Community Standard (D6300).

1 Vortex Genie with 2ml BashingBead™ Tubes

Recommended for ease of use and accessibility

Use Microtube Adaptor (Scientific Industries, Inc. Cat. No. S5001-7)

1. 40 minutes of continuous bead beating (max of 18 tubes per adaptor)

2 Bertin Precellys Evolution with 2 ml BashingBead™ Tubes

Recommended for ease of use and ultra-high speed.

1. 1 minute on at 9,000 RPM
2. 2 minutes rest
3. Repeat cycle 4 times for a total of 4 minutes of bead beating

3 MP Fastprep-24™ (Classic & 5G) with 2 ml BashingBead™ Tubes

Maximum of 20 tubes. The weight of > 20 tubes may cause a system error.

1. 1 minute on at 6.5 m/s
2. 5 minutes rest
3. Repeat cycle 5 times for a total of 5 minutes of bead beating

4 Omni Bead Ruptor Elite with 2 ml BashingBead™ Tubes

1. 1 minute on at 6 m/s
2. 5 minutes rest
3. Repeat cycle 3 times for a total of 3 minutes of bead beating

5 Biospec Mini-BeadBeater-16 with 2 ml BashingBead™ Tubes

1. 1 minute at maximum speed
2. 5 minutes rest
3. Repeat cycle 5 times for a total of 5 minutes of bead beating

6 Biospec Mini-BeadBeater-96 with 2 ml BashingBead™ Tubes

1. 5 minutes on at Max RPM
2. 5 minutes rest
3. Repeat cycle 4 times for a total of 20 minutes of bead beating

7 Biospec Mini-BeadBeater-96 with 96 well lysis rack

1. 5 minutes on at Max RPM
2. 5 minutes rest
3. Repeat cycle 8 times for a total of 40 minutes of bead beating

✗ TissueLyser II

No tested conditions yielded accurate profiles. This device is not validated by Zymo Research for microbiome research.

✗ TissueLyser LT

No tested conditions yielded accurate profiles. This device is not validated by Zymo Research for microbiome research.

✗ Retsch Mixer Mill MM 400

No tested conditions yielded accurate profiles. This device is not validated by Zymo Research for microbiome research.