Fecal Collection Kit

WARNING: DO NOT drink or touch the solution. The solution can be harmful if ingested and may irritate the skin. Avoid contact with eyes. For specific information, including instructions on how to use the kit, please refer to the product safety Data Sheet (SDS).

15°C / 25°C

DNA/RNA Shield Fecal Collection Kit

REF R1180
For At-Home Sample Collection

English User Instructions

Kit Contents:
1 x Fecal Tube
1 x FecesCatcher
1 x Biohazard Bag
1 x Gloves
1 x Instructions

READ ALL INSTRUCTIONS PRIOR TO USE

Summary & Explanation of the Kit

This kit provides the instructions and materials for collecting and stabilizing fecal specimens.

Intended Use

For the collection of fecal samples for DNA and RNA applications.

For research use only
Not for use in diagnostic applications

Contents

Kit contains tube pre-filled with stabilizing solution, gloves, FecesCatcher, biohazard bag, and instructions for use.

Collection Precaution

Do NOT spill or ingest the stabilizing solution inside the tube. Wash with water if stabilizing solution comes in contact with eyes or skin.

Procedure

Note: Prevent feces from falling into toilet water to avoid sample contamination.

1. Open the FecesCatcher carefully and follow instructions printed on the catcher for placing onto toilet seat.
2. After bowel movement, unscrew the tube cap. Be careful not to spill or ingest solution inside.
3. Use the spoon to scoop one spoonful of feces (approximately the size of a pea).
4. Place the collected fecal sample into the collection tube and screw tightly.
5. Ensure cap is tight and shake to mix for 10 seconds.
6. After collection, detach the FecesCatcher from toilet seat and flush.

Storage: 15°C / 25°C

WARNING

FOR EXTERNAL USE ONLY. DO NOT drink, touch or remove the solution from the tube. Avoid skin and eye contact with the liquid in the tube. DO NOT ingest. IF SWALLOWED: Call a doctor if you feel unwell. Rinse mouth. IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. SMALL CAP MAY POSE A CHOKING HAZARD.

Label Legend

- Collect Sample By
- Catalog Number
- Manufacturer
- Storage Instructions
- Harmful
- Lot Number
- See Instruction Manual

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
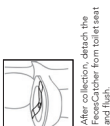
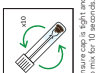
1. Open the FecesCatcher carefully and follow instructions printed on the catcher for placing onto toilet seat.

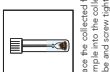
2. After-bowel movement, unscrew the tube cap. Be careful not to spill or ingest solution inside.

3. Use the spoon to scoop one spoonful of feces (approximately the size of a pea).





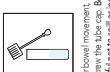
4. Place the collected fecal sample into the collection tube and screw tightly.

5. Ensure cap is tight and shake to mix for 10 seconds.

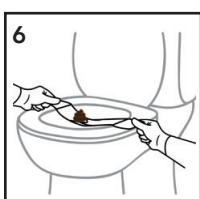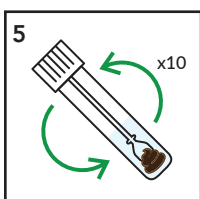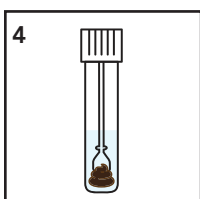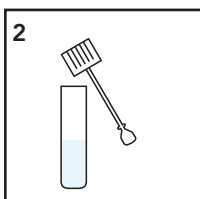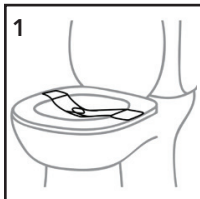6. After collection, detach the FecesCatcher from toilet seat and flush.

Scan the QR code for video instructions

Wash with water if stabilization reagent comes in contact with eyes or skin. DO NOT ingest.